

Human IFN γ ELISpot Matched Antibody Pair

Catalog No. CDK208

Quantity: 10 x 96 wells

Intended Use

The Cell Sciences® ELISpot is a highly specific immunoassay for the analysis of cytokine and other soluble molecule production and secretion from T-cells at a single cell level in conditions closely comparable to the in-vivo environment with minimal cell manipulation. This technique is designed to determine the frequency of cytokine producing cells under a given stimulation and the comparison of such frequency against a specific treatment or pathological state. The ELISpot assay constitutes an ideal tool in the investigation of Th1 / Th2 responses, vaccine development, viral infection monitoring and treatment, cancerology, infectious disease, autoimmune diseases and transplantation.

Utilizing sandwich immuno-enzyme technology, Cell Sciences® ELISpot assays can detect both secreted cytokines and single cells that simultaneously produce multiple cytokines. Cell secreted cytokines or soluble molecules are captured by coated antibodies avoiding diffusion in supernatant, protease degradation or binding on soluble membrane receptors. After cell removal, the captured cytokines are revealed by tracer antibodies and appropriate conjugates.

This kit has been configured for research use only and is not to be used in diagnostic procedures.

2. Introduction

2.1. Summary

Different populations of T-cells secrete differing patterns of cytokines that ultimately lead to different immune responses. IFN γ production is a key function of Th1, CD8⁺ CTLs and also NK cells. IFN γ is a cytokine critical for cell mediated immunity against viral and intracellular bacterial infections and is involved in the inflammatory response following secretion via macrophage activation and stimulation of antibody secretion. IFN γ is the hallmark effector cytokine of Th1 and therefore is an excellent marker for identifying a host response to intracellular pathogens.

IFN γ is produced during infection by T cells of the cytotoxic/suppressor phenotype (CD8) and by a subtype of helper T cells, the Th1 cells. Th1 cells secrete IL-2, IL-3, TNF α and IFN γ , whereas Th2 cells mainly produce IL-3, IL-4, IL-5, and IL-10, but little or no IFN γ . IFN- γ preferentially inhibits the proliferation of Th2 but not Th1 cells, indicating that the presence of IFN γ during an immune response will result in the preferential proliferation of Th1 cells.

In addition, IFN- γ has several properties related to immunoregulation. IFN γ is a potent activator of mononuclear phagocytes, and activates macrophages to kill tumor cells by releasing reactive oxygen intermediates and TNF α . IFN γ induces or augments the expression of MHC antigens on macrophages, T and B cells and some tumor cell lines. On T and B cells IFN- γ promotes differentiation. It enhances proliferation of activated B cells and can act synergistically with IL-2 to increase immunoglobulin light-chain synthesis.



Cell Sciences, Inc.
480 Neponset Street
Bldg 12A
Canton, MA 02021


Toll Free: 888-769-1246
Phone: 781-828-0610
Fax: 781-828-0542

E-mail: info@cellsciences.com
Web Site: www.cellsciences.com

The role of IFN γ as a disease marker has been demonstrated for a number of different pathological situations including, viral infection, Autoimmune disease, transplant rejection, Diabetes and allergy.

2.2. Principle of the method

Capture antibodies highly specific for the analyte of interest are coated to the wells of a PVDF bottomed 96 well microtiter plate either during kit manufacture or in the laboratory. The plate is then blocked to minimize any non-antibody dependent unspecific binding and finally washed before adding the cells to be investigated. Cell suspension and stimulant are added to the coated and blocked microtiter plate and the plate incubated allowing the specific antibodies to bind any analytes produced. Biotinylated detection antibodies are then added which bind to the previously captured analyte. Enzyme conjugated streptavidin is added binding to the detection antibodies. Any excess unbound analyte and antibodies are removed by careful washing. Color substrate is then applied to the wells resulting in colored spots, which can be quantified using appropriate analysis software or manually using microscopes.

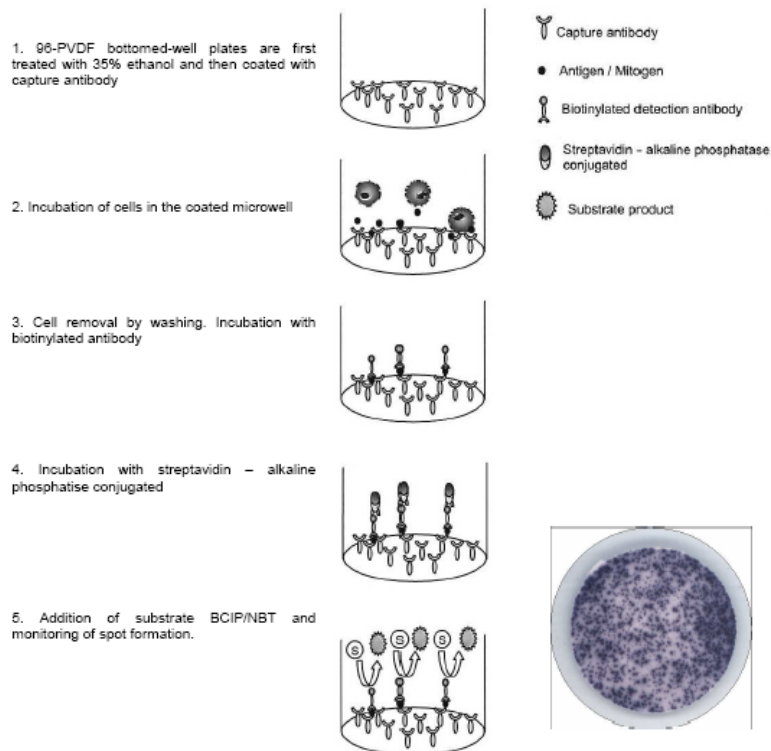

3. Reagents provided (for 10 x 96 tests)

- Capture Antibody (2 vials of 0.5ml). The antibody is supplied sterile and does not contain preservative. We strongly advise sterile pipetting.
- Biotinylated detection antibody (2 vials, lyophilized)



4. Materials/Reagents required but not provided

- Miscellaneous laboratory plastic and/or glass, if possible sterile
- Streptavidin-Alkaline Phosphates conjugated
- Bovine Serum Albumin (BSA)
- Substrate solution (BCIP/NBT)
- Ethanol
- Cell culture reagents (e.g. RPMI-1640, L-glutamine, and FCS)
- Cell stimulation reagents (e.g. PMA and Ionomycin)
- CO2 incubator
- Tween 20
- Phosphate Buffered Saline (PBS)
- 96 well PVDF bottomed plates

5. Storage Instructions

Store kit reagents between 2 and 8°C except uncoated plates, which should be stored at RT. Immediately after use, remaining reagents should be returned to cold storage (2 to 8°C). Expiry of the kit and reagents is stated on box front labels. The expiry of the kit components can only be guaranteed if the components are stored properly, and if in the case of repeated use of one component, the reagent is not contaminated by the first handling.

6. Safety & Precautions for use

- For research use only not to be used as a diagnostic test
- Handling of reagents, serum or plasma specimens should be in accordance with local safety procedures, e.g. CDC /NIH Health manual: "Biosafety in Microbiological and Biomedical Laboratories" 1984
- Do not eat, drink, smoke or apply cosmetics where kit reagents are used
- Do not pipette by mouth
- When not in use, kit components should be stored refrigerated or frozen as indicated on vials or bottles labels
- All reagents should be warmed to room temperature before use.
- Cover or cap all reagents when not in use
- Do not mix or interchange reagents between different lots
- Do not use reagents beyond the expiration date of the kit
- Use a clean disposable plastic pipette tip for each reagent, standard, or specimen addition in order to avoid cross contamination
- Use a clean plastic container to prepare the washing solution
- Thoroughly mix the reagents and samples before use by agitation or swirling
- All residual washing liquid must be drained from the wells by efficient aspiration or by decantation followed by tapping the plate forcefully on absorbent paper. Never insert absorbent paper directly into the wells
- When pipetting reagents, maintain a consistent order of addition from well-to-well. This will ensure equal incubation times for all wells
- **BCIP/NBT buffer** is potentially carcinogenic and should be disposed of appropriately, caution should be taken when handling this reagent, always wear gloves
- Follow incubation times described in the assay procedure



7. Reagent Preparation

7.1. 1X Phosphate Buffered Saline (PBS) (Coating Buffer)

For 1 liter of 10X PBS weigh-out: 80g NaCl
2g KH₂PO₄
14.4g Na₂HPO₄ · 2H₂O.

Add distilled water to 1 liter. Adjust the pH of the solution to 7.4 +/- 0.1 were required.

Dilute the solution to 1X before use.

7.2. Cell culture media + 10% Serum (Blocking Buffer)

For one plate add 1ml Serum (e.g. FCS) to 9ml of culture media (use same cell culture medium as used to derive the cell suspension).

7.3. 1% BSA PBS Solution (Dilution Buffer)

For one plate, dissolve 0.2 g of BSA in 20 ml of 1X PBS.

7.4. 0.05% PBS-T Solution (Wash Buffer)

For one plate, dissolve 50µl of Tween 20 in 100 ml of 1X PBS.

7.5. 35% Ethanol (PVDF Membrane Activation Buffer)

For one plate, mix 3.5 ml of ethanol with 6.5 ml of distilled water.

7.6. Capture Antibody

This reagent is supplied sterile once opened keep the vial sterile or aliquot and store at -20oC. For optimal performance prepare the Capture Antibody dilution immediately before use.

Dilute 100µl of capture antibody in 10 mL of 1X PBS and mix well.

7.7. Detection Antibody

Reconstitute the lyophilized antibody with 0.55mL of distilled water. Gently mix the solution and wait until all the lyophilized material is back into solution.

If not used within a short period of time, reconstituted Detection Antibody should be aliquoted and stored at -20°C. In these conditions the reagent is stable for r at least one year. For optimal performance prepare the reconstituted antibody dilution immediately prior to use.

Dilute 100ml of antibody into 10ml Dilution Buffer and mix well.

7.8. Streptavidin – Alkaline Phosphates

Dilute in Dilution buffer. DO NOT KEEP THE DILUTIONS FOR FURTHER EXPERIMENTS.



8. Sample and Control Preparation

8.1. Cell Stimulation

Cells either can be stimulated directly in the antibody coated wells (Direct) or, first stimulated in 24 well plates or flask, harvested, and then plated into the coated wells (Indirect).

The method used is dependent on 1) the type of cell assayed 2) the expected cell frequency. When a low number of cytokine producing cells are expected it is also advised to test them with the direct method, however, when this number is particularly high it is better to use the indirect ELISpot method.

All the method steps following stimulation of the cells are the same whatever the method (direct/indirect) chosen.

8.2. Positive Assay Control, IL-10 production

We recommend using the following polyclonal activation as a positive control in your assay.

Dilute PBMC in culture media (e.g. RPMI 1640 supplemented with 2mM L-glutamine and 10% heat inactivated fetal calf serum) containing 1ng/ml PMA and 500ng/ml Ionomycin (Sigma, Saint Louis, MO). Distribute 2.5×10^4 to 5×10^4 cells per 100 μ l in required wells of an antibody coated 96-well PVDF plates and incubate for 15-20 hours in an incubator.

For other stimulators incubation times may vary, depending on the frequency of cytokine producing cells, and should be optimized in each situation.

8.3. Negative Assay Control

Dilute PBMC in culture media to give an appropriate cell number (same number of un-stimulated cells as stimulated sample cells) per 100 μ l with no stimulation.

8.4. Sample

Dilute PBMC in culture medium and stimulator of interest (i.e. Sample, Vaccine, Peptide pool or infected cells) to give an appropriate cell number per 100 μ l.

Optimal assay performances are observed between 1×10^5 and 2.5×10^5 cells per 100ml.

Stimulators and incubation times can be varied depending on the frequency of cytokine producing cells and therefore should be optimized by the testing laboratory.



9. Method

Prepare all reagents as shown in section 7 and 8.

Note: For optimal performance, prepare the Streptavidin-AP dilution immediately prior to use

Assay Step	Details
1. Addition	Add 25µl of 35% ethanol to every well
2. Incubation	Incubate plate at room temperature (RT) for 30 seconds
3. Wash	Empty the wells by flicking the plate over a sink & gently tapping on absorbent paper. Thoroughly wash the plate 3x with 100µl of 1X PBS per well
4. Addition	Add 100µl of diluted capture antibody to every well
5. Incubation	Cover the plate and incubate at 4°C overnight
6. Wash	Empty the wells as previous and wash the plate once with 100µl of 1X PBS per well
7. Addition	Add 100µl of culture media with 10% serum to every well
8. Incubation	Cover the plate and incubate at RT for 2 hours
9. Wash	Empty the wells as previous and thoroughly wash once with 100µl of 1X PBS per well
10. Addition	Add 100µl of sample, positive and negative controls cell suspension to appropriate wells providing the required concentration of cells and stimulant (cells may have been previously stimulated see section 8.)
11. Incubation	Cover the plate and incubate at 37°C in a CO ₂ incubator for an appropriate length of time (15-20 hours). Note: do not agitate or move the plate during this incubation
12. Addition	Empty the wells and remove excess solution then add 100µl of PBS-T to every well
13. Incubation	Incubate the plate at 4°C for 10 min
14. Wash	Empty the wells as previous and wash the plate 3x with 100µl of PBS-T
15. Addition	Add 100µl of diluted detection antibody to every well
16. Incubation	Cover the plate and incubate at RT for 1 hour 30 min
17. wash	Empty the wells as previous and wash the plate 3x with 100µl of PBS-T
18. Addition	Add 100µl of diluted Streptavidin-AP conjugate to every well
19. Incubation	Cover the plate and incubate at RT for 1 hour
20. Wash	Empty the wells and wash the plate 3x with 100µl of PBS-T
21. Wash	Peel of the plate bottom and wash both sides of the membrane 3x under running distilled water, once washing complete remove any excess solution by repeated tapping on absorbent paper.
22. Addition	Add 100µl of ready-to-use BCIP/NBT buffer to every well
23. Development	Incubate the plate for 5-15 min monitoring spot formation visually throughout the incubation period to assess sufficient colour development
24. Wash	Empty the wells and rinse both sides of the membrane 3x under running distilled water. Completely remove any excess solution by gentle repeated tapping on absorbent paper
<p>Read Spots: allow the wells to dry and then read results. The frequency of the resulting coloured spots corresponding to the cytokine producing cells can be determined using an appropriate ELISpot reader and analysis software or manually using a microscope.</p> <p><i>Note: spots may become sharper after overnight incubation at 4°C</i></p>	

Plate should be stored at RT away from direct light, but please note color may fade over prolonged periods so read results within 24 hours.



10. Performance Characteristics

10.1. Specificity

The assay recognizes natural human IFN γ .

To define specificity of this ELISpot, several proteins were tested for cross reactivity. There was no cross reactivity observed for these proteins tested (IL-1 α , IL-1 β , IL-10, IL-12, TNF α , IL-8, and IL-13. This testing was performed using the equivalent IFN γ antibody pair in an ELISA assay.

NOT FOR HUMAN USE. FOR RESEARCH ONLY. NOT FOR DIAGNOSTIC OR THERAPEUTIC USE.



Cell Sciences, Inc.
480 Neponset Street
Bldg 12A
Canton, MA 02021

Toll Free: 888-769-1246
Phone: 781-828-0610
Fax: 781-828-0542

E-mail: info@cellsciences.com
Web Site: www.cellsciences.com